


User Guide

Mantis Elite-Cam HD Connection & Setup

Mantis Elite-Cam HD Integrated USB Digital Camera

Mantis Elite-Cam HD is a permanently-engaged USB digital camera solution for use in a wide range of tasks requiring magnification.

Health & Safety

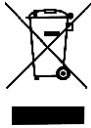

Vision Engineering and its products conform to the requirements of the EC Directives on Waste Electrical and Electronic Equipment (WEEE) and Restriction of Hazardous Substances (RoHS).

INTRODUCTION

Mantis Elite-Cam HD Integrated USB Digital Camera	i
Health & Safety	i

CONTENTS

Pack contents	1
---------------	---


GETTING STARTED

Assembly	2
Operation	2
General care and maintenance	2

WARRANTY


Pack contents

- 1 Mantis Elite-Cam HD viewing head
- 2 Software USB memory stick
- 3 Interconnect cable
- 4 Mantis Elite-Cam HD software user guide (not shown)
- 5 Mantis Elite user guide (not shown)
- 6 Mantis Elite-Cam HD user guide (not shown)

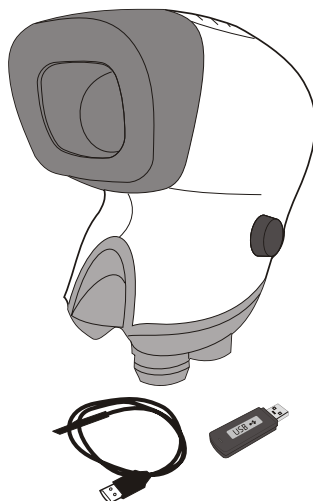

Assembly

- ▶ Follow the Mantis Elite user guide to mount the viewing head **1** to the stand.
- ▶ Install software contained on the software USB stick **2**, as per Mantis Elite-Cam HD software user guide.
- ▶ Before connecting to the chosen computer install the uEye software as detailed in the Mantis Elite-Cam HD software user guide.
- ▶ Connect cable **3** to the viewing head **1** and then the chosen computer.

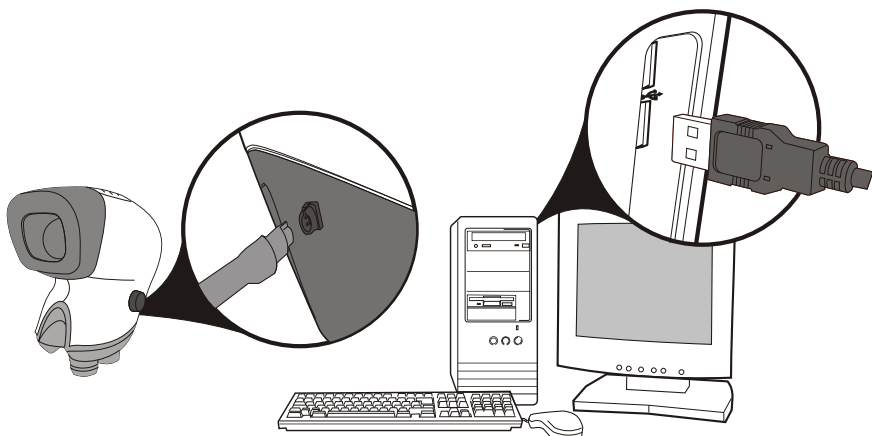

Operation

- ▶ Refer to Mantis Elite user guide provided for general operation.
- ▶ Refer to the Mantis Elite-Cam HD software user guide for details on how to start using the supplied uEye software.

General care and maintenance

For all care and maintenance guidelines see the Mantis Elite user guide.



WARRANTY

This product is warranted to be free from defects in material and workmanship for a period of one year from the date of invoice to the original purchaser.

If during the warranty period the product is found to be defective, it will be repaired or replaced at facilities of Vision Engineering or elsewhere, all at the option of Vision Engineering. However, Vision Engineering reserves the right to refund the purchase price if it is unable to provide replacement, and repair is not commercially practicable or cannot be timely made. Parts not of Vision Engineering manufacture carry only the warranty of their manufacturer. Expendable components such as fuses carry no warranty.

This warranty does not cover damage in transit, damage caused by misuse, neglect, or carelessness, or damage resulting from either improper servicing or modification by other than Vision Engineering approved service personnel. Further, this warranty does not cover any routine maintenance work on the product described in the user guide or any minor maintenance work which is reasonably expected to be performed by the purchaser.

No responsibility is assumed for unsatisfactory operating performance due to environmental conditions such as humidity, dust, corrosive chemicals, deposition of oil or other foreign matter, spillage, or other conditions beyond the control of Vision Engineering.

Except as stated herein, Vision Engineering makes no other warranties, express or implied by law, whether for resale, fitness for a particular purpose or otherwise. Further, Vision Engineering shall not under any circumstances be liable for incidental, consequential or other damages.

For more information...

Vision Engineering has a network of offices and technical distributors around the world. For more information, please contact your Vision Engineering branch, local authorised distributor, or visit our website.

Distributor



Vision Engineering Ltd.
(Manufacturing)
Send Road, Send,
Woking, Surrey GU22 7ER, England
Tel: +44 (0) 1483 248300
Email: generalinfo@visioneng.com

Vision Engineering Ltd.
(Commercial)
Monument House, Monument Way West,
Woking, Surrey, GU21 5EN, England
Tel: +44 (0) 1483 248300
Email: generalinfo@visioneng.com

Vision Engineering Inc.
(Manufacturing & Commercial)
570 Danbury Road,
New Milford, CT 06776, USA
Tel: +1 (860) 353 3776
Email: info@visioneng.com

Vision Engineering Inc.
(West Coast Commercial)
745 West 34th Avenue,
Cheng, CA 92605, USA
Tel: +1 (714) 974 6966
Email: info@visioneng.com

Vision Engineering Ltd.
(Central Europe)
Anton-Pendele-Str 3,
82275 Emming, Deutschland
Tel: +49 (0) 8141 40167-0
Email: info@visioneng.de

Vision Engineering Ltd.
(France)
ZAC de la Némésais, Av. de la Némésais
91220 Le Plessis Palé, France
Tel: +33 (0) 160 76 40 00
Email: info@visioneng.fr

Vision Engineering Ltd.
(Italy)
Via Cesare Canto, 9
20092 Cinisello Balsamo MI, Italia
Tel: +39 (0) 4729 3518
Email: info@visioneng.it

Vision Engineering
(Brazil)
Email: info@visioneng.com.br

Nippon Vision Engineering
(Japan)
272-2 Saedo-cho, Tsudaki-ku,
Yokohama-shi, 224-0054, Japan
Tel: +81 (0) 45 935 1117
Email: info@visioneng.jp

Vision Engineering
(China)
111, International Ocean Building,
720 Pudong Avenue, Shanghai,
200120, P.R. China
Tel: +86 (0) 21 5036 7556
Email: info@visioneng.com.cn

Vision Engineering
(S.E. Asia)
Tel: +603 80700908
Email: info@visioneng.asia

Vision Engineering
(India)
Email: info@visioneng.co.in

Visit our website:

www.visioneng.com

Disclaimer – Vision Engineering Ltd. has a policy of continuous development and reserves the right to change or update, without notice, the design, material or specification of any products. The information contained within this brochure/datasheet and to discontinue production or distribution of any of the products described.